

# Direct “on-resin” radiofluorination, radiotitanation and radiozirconation using automated radiosynthesis module BioSyntheSizer 2.1

Fedor Zhuravlev†\*, David Tezét†, Jesper Fonslett†, Toni Bohatzsch§, and Steffen Howitz§



†Hevesy Laboratory, Center for Nuclear Technologies at the Technical University of Denmark, Frederiksborgvej 399, 4000 Roskilde, Denmark.

§GeSiM mbH, Bautzner Landstrasse 45, D-01454 Radeberg, Germany

\*Correspondence email: fezb@dtu.dk

## Introduction

Although Nuclear Medicine market is dominated by Diagnostics, Research segment grows by an impressive 10.2% CAGR. ▶

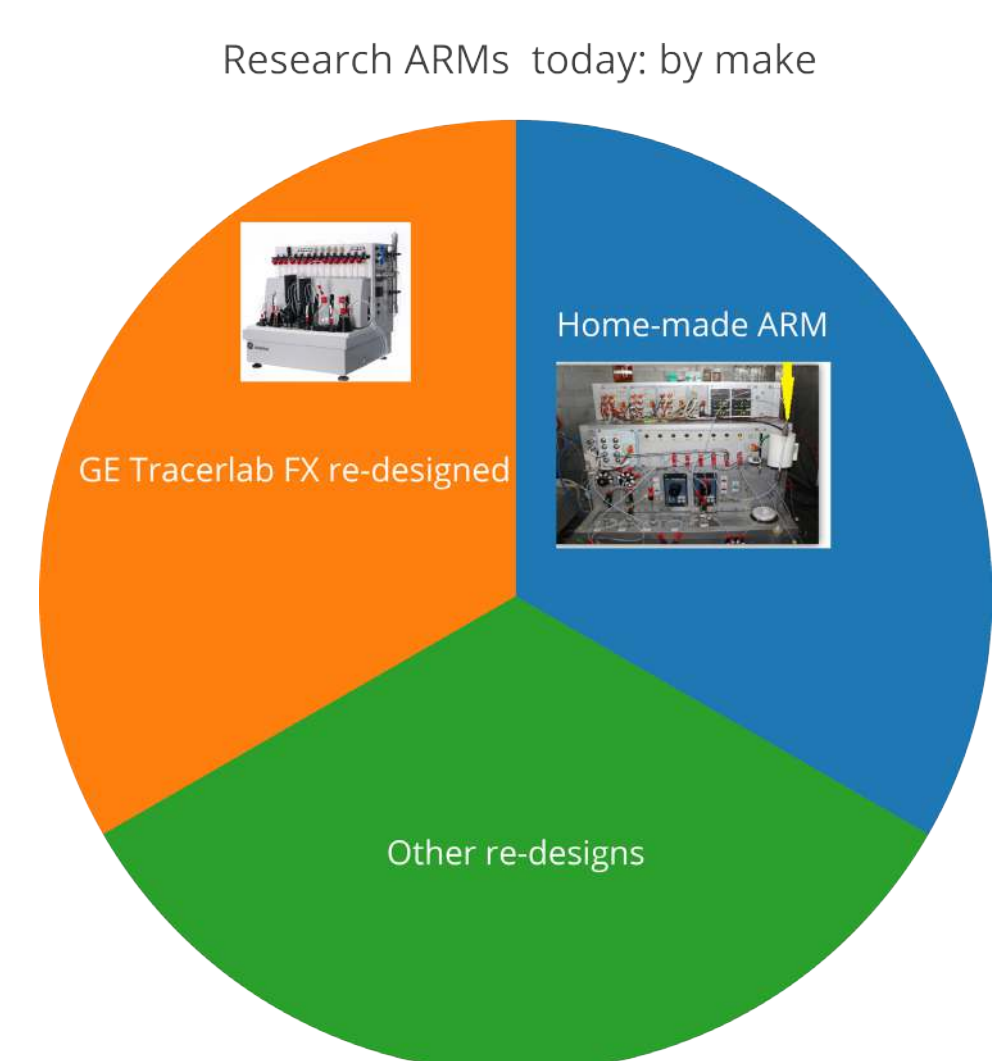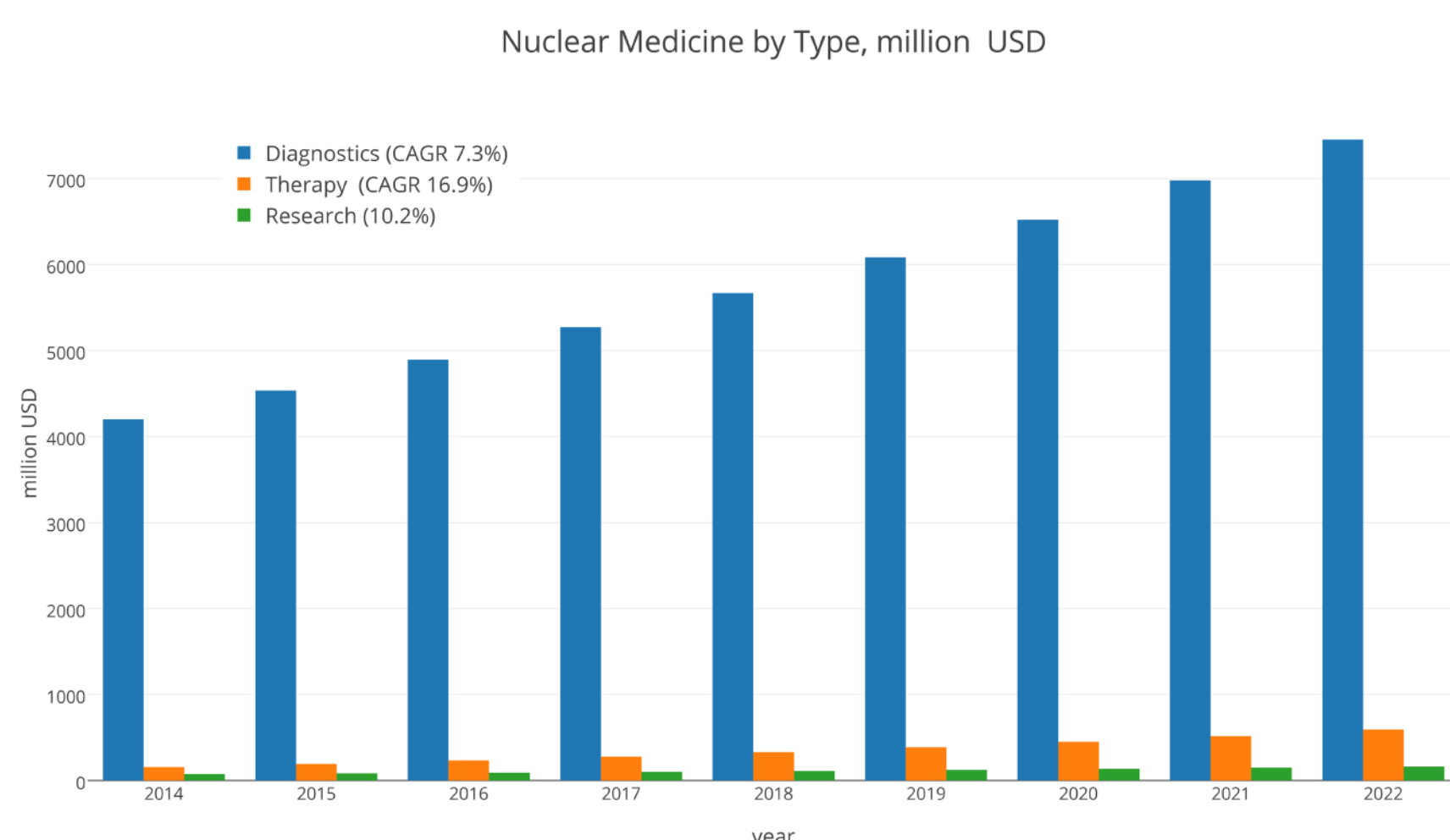

There is an unmet need for automated radiosynthesis modules (ARMs), catering specifically to R&D segment of the market.

◀ According to our survey, home-made or re-designed ARMs are currently the most popular solutions for R&D labs.

## Objectives

Our research program is directed towards new approaches to automated radiofluorination and radiometallation using radiolabeling at the point of radionuclide capture [1-2]. Herein we report automated trapping, trap-and-release, and “on-resin” radiolabeling with  $^{18}\text{F}$ ,  $^{45}\text{Ti}$ , and  $^{89}\text{Zr}$  using radiosynthesis module BioSyntheSizer 2.1 (GeSiM mbH, Germany).

## BioSyntheSizer platform

BioSyntheSizer 2.1 (BSys) is an advanced robotic platform for complex laboratory automation. A toolhead movable in 3D controls fluidics via luer-lock manifolds connected to syringe pumps for easy aspiration/dispensing operations. The reagent rack can host a variety of vials (1-50 mL), and solid-phase extraction (SPE) cartridges.

A unique feature is handling of syringe needles and Eppendorf pipette tips.



◀ A 3D movable toolhead combines the versatility of manual operator with reproducibility of a machine.


## Solid Phase Flow Reaction Module

Our developer’s version of BSys is equipped with a solid phase flow reaction module. The module is electronically integrated and software-controlled in the BSys allowing for trapping, trap-and-release, and “on-resin” radiolabeling.

The module features:

1. Fine temperature control
2. Accommodates different sizes of glass columns for radiofluorination and radiometallation

## Radiofluorination

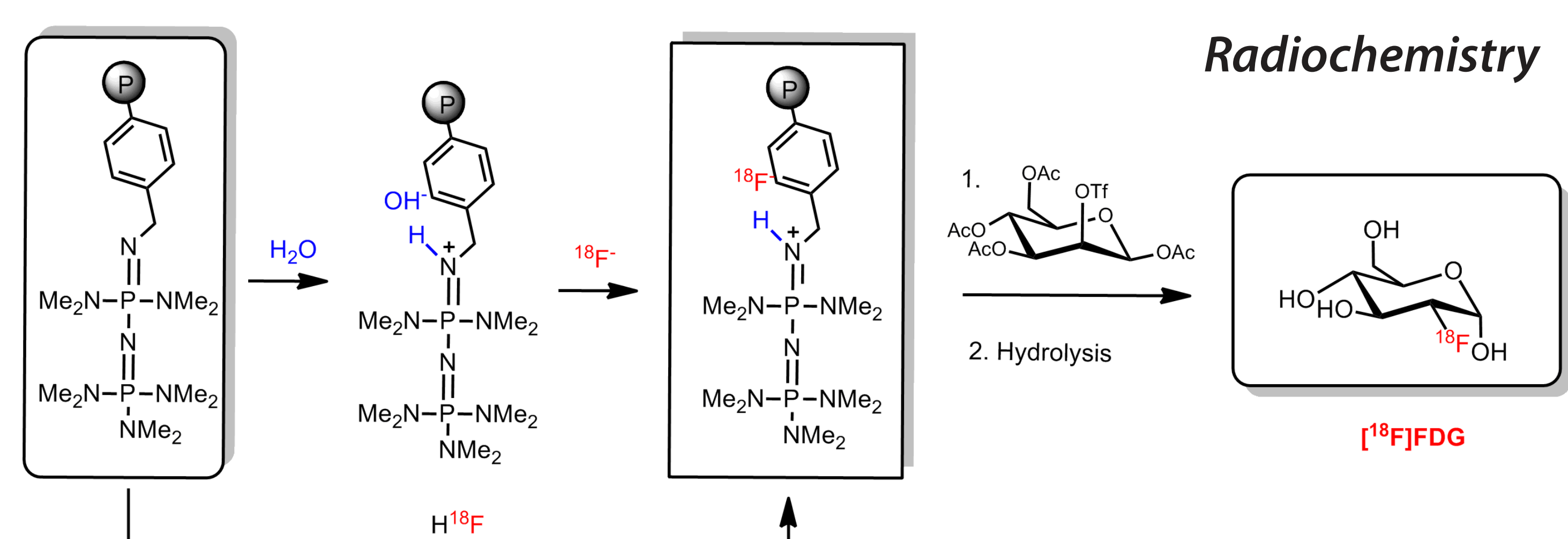

1. Column is loaded with polymer-supported phosphazene
2. Target water is passed through, (L1→L2) 98-99% of fluoride is trapped
3. A quick wash with acetonitrile (ACN) removes residual water from the column (does not have to be very dry!)
4. A substrate in a suitable solvent (toluene, ACN) is passed through the column while the column is heated at 80 - 100 °C (L1→L2)
5. The product is collected in L2. RCY = 40–60%.
6. With many substrates the resin can be reused at least 4 times

## Radiometallation: $^{45}\text{Ti}$ , $^{89}\text{Zr}$

$^{45}\text{Ti}$  is great for peptide & affibody PET


$^{89}\text{Zr}$  is great for immuno PET

- 3.08 h half life
- 85% positron branch
- $E_0 = 1.04 \text{ MeV}$
- $E_{\text{ave}} = 439 \text{ keV}$
- $^{nat}\text{Sc}(p,n)^{45}\text{Ti}$

- 78.4 h half life
- 23% positron branch
- $E_0 = 0.90 \text{ MeV}$
- $E_{\text{ave}} = 396 \text{ keV}$
- $^{nat}\text{Y}(p,n)^{89}\text{Zr}$



## Targetry

23 mg Sc  
8 Gbq EOB

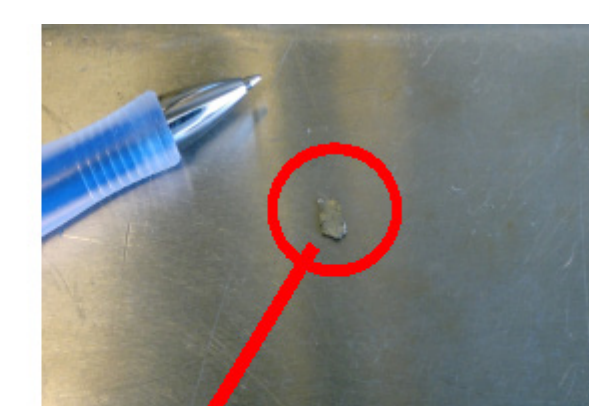


## Automated target foil digestion: no manual handling

(A)  $^{45}\text{Ti}$ ,  $^{89}\text{Zr}$ : 37% HCl  
(B)  $^{89}\text{Zr}$ : 4M HCl/dioxane iPrOH

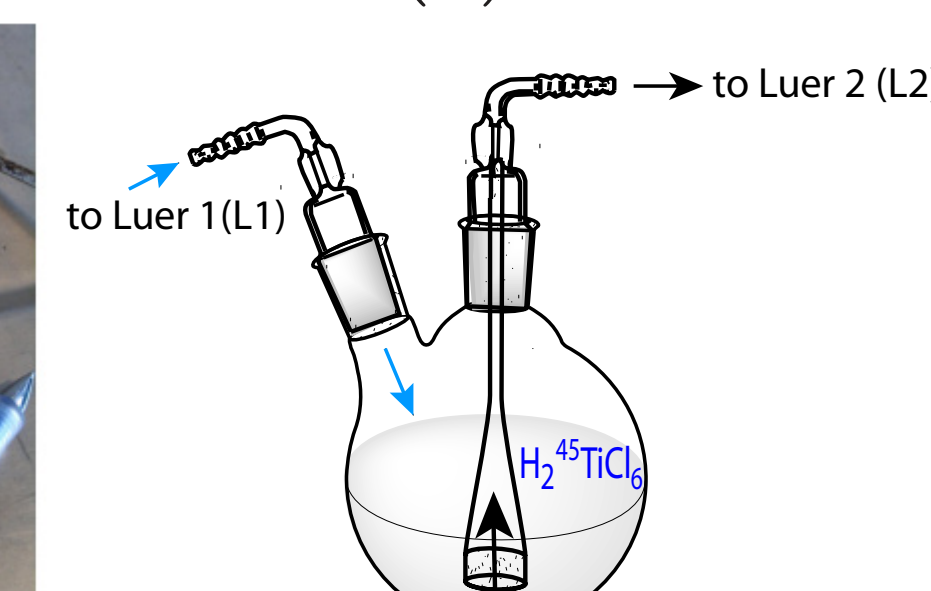
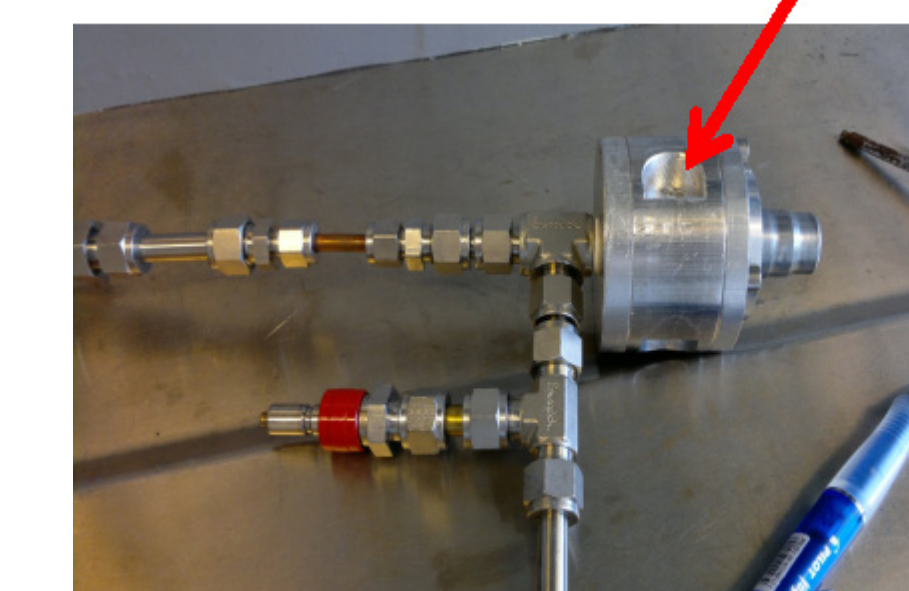


Automated digestion: ▶ [YouTube](#)

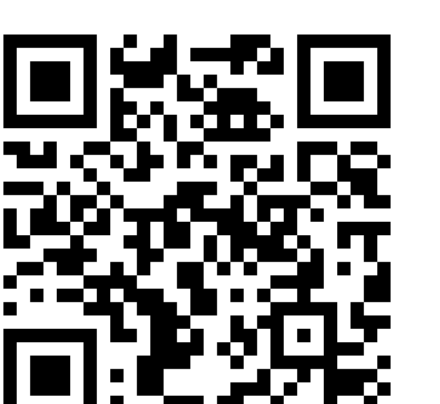

## Radiochemistry

Distribution constants:  $K^D_{\text{Ti}} = 91.4, \pm 10.5 \text{ mL/mg}$  and  $K^D_{\text{Sc}} = 6.8 \pm 1.4 \text{ mL/mg}$

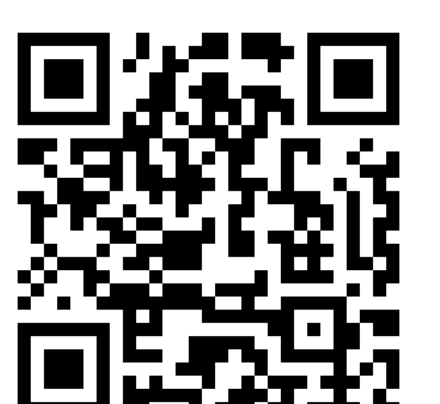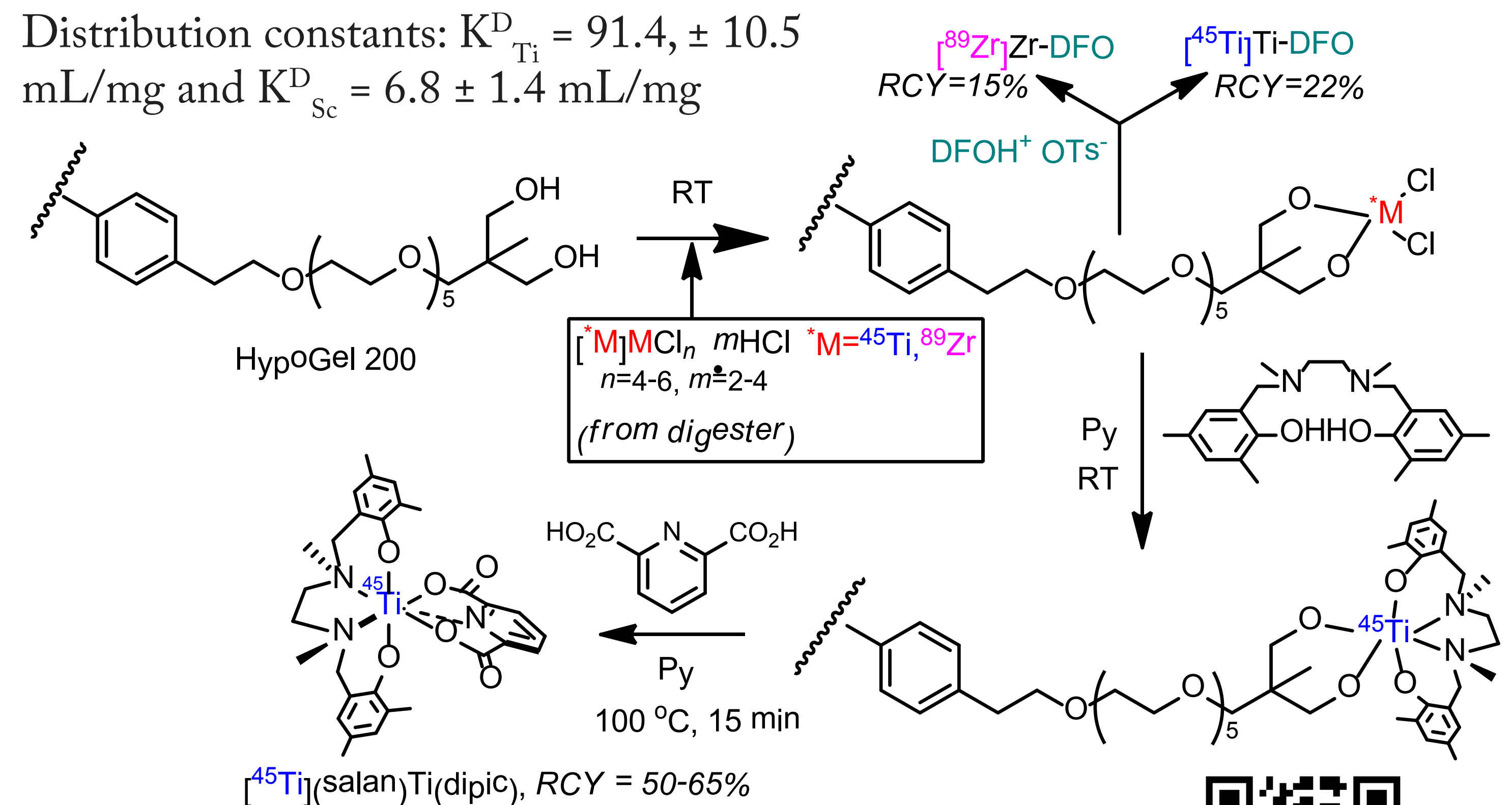

## Conclusions

1. Integrated target digester and the solid phase flow reactor enabled automated radiosynthesis of  $^{18}\text{F}$ -FDG, and showcased radiometallation with  $^{45}\text{Ti}$  and  $^{89}\text{Zr}$  directly on a radionuclide trapping resin.
2. The fluidics and robotics proved to be tolerant to fuming HCl, and non-aqueous HCl/dioxane.
3. The flexible design of the radiosynthesizer allowing for creation of fluid path on demand via a 3D movable fluidics head required no hardware re-configuration in-between radiometal and fluorine-18 radiolabeling

**Acknowledgements:** David Tezét was supported by the Proof of Concept grant from DTU, Denmark.

**References:** [1] Mathiessen, B.; Zhuravlev, F.\* *Molecules* **2013**, *18* (9), 10531–10547.

[2] Severin, G. W.; Nielsen, C. H.; Jensen, A. I.; Fonslet, J.; Kjær, A.; Zhuravlev, F.\* *J. Med. Chem.* **2015**, *58* (18), 7591–7595.